



Poster Reprint

Controlled Release Society

July 2007

The Influence of Plasticizer Type and Concentration on Acid Resistance of Tablets Coated with a New Aqueous Delayed Release Film Coating System

Cara Young, Scott Vass, Kurt Fegely, and Ali Rajabi-Siahboomi
Colorcon, Inc., Moyer Boulevard, West Point, Pennsylvania, USA,

Abstract Summary

The influence of plasticizer type and concentration, added to a new aqueous delayed-release (DR) film coating system, on acid resistance was evaluated. Incorporating PEG 8000 (8-15%), TEC (12-25%), or triacetin (12-30%) in the DR coating system resulted in acid resistance (2 hours) and complete disintegration in buffer (pH 6.8), as desired.

Introduction

Enteric protection from DR film coating systems has been shown to be influenced by the hydrophilic and lipophilic properties of the plasticizer, level of plasticizer, and amount of polymer applied to the substrate¹. Furthermore, the properties of the substrate, in conjunction with the plasticizer type, were found to influence the degree of enteric protection².

Therefore, the purpose of this study was to evaluate the influence of three plasticizers at various levels on acid resistance in a new aqueous DR film coating system (Acryl-EZE[®] 93A Series). The tablets were tested in two different media, 0.1N HCl (pH 1.2) and 0.5M sodium acetate buffer (pH 4.5) to simulate the stomach pH of a fasted human and the increase in stomach pH (above pH 4.0) from food and/or multiple doses of proton pump inhibitors (PPIs), respectively. A multiple dose regimen of PPIs results in a decrease in gastric acid secretion with a subsequent elevation in gastric PH⁴⁻⁵.

The properties of the substrate (placebo cores) and amount of polymer applied to the substrate (7.6mg/cm²) were held constant. Polyethylene glycol 8000 (8-15%), triethyl citrate (12-25%), or triacetin (12-30%) was added to water, followed by Acryl-EZE 93A, to form a pigmented aqueous DR coating dispersion. Placebo tablets were coated with the DR coating dispersion to 10% theoretical weight gain. Acid uptake in 0.1N HCl and acetate buffer (pH 4.5), as well as disintegration time in phosphate buffer (pH 6.8), were determined.

Experimental Methods

Materials

Polyethylene glycol 8000 (PEG 8000), triethyl citrate (TEC), and triacetin were obtained from Clariant (USA), Morflex (USA), and Tessenderlo Fine Chemicals (UK), respectively. Simethicone emulsion (30% Dow) was used as an anti-foam to de-aerate the DR coating dispersion, was purchased from Dow (USA). Placebo tablets, Opadry[®] YS-1-7027, and Acryl-EZE 93A18597 were obtained from Colorcon (USA).

Preparation and Characterization

A seal coat dispersion was prepared by dispersing (low shear) Opadry in water and mixing for 30 minutes. The solids content of the final coating dispersion was 12.5% (w/w).

The enteric coating dispersion was prepared by dispersing the plasticizer and simethicone emulsion (0.5% w/w, respect to total dispersion) in water and mixing (low shear) for 5 minutes. Acryl-EZE 93A was added (20% solids, w/w, respect to water) and mixed for an additional 30 minutes, achieving a uniform DR pigmented coating dispersion. The dispersion was screened (250µm) and slowly stirred during spraying.

Placebo tablets (3/8" (9.57mm) convex, 307mg) were seal coated (4% theoretical weight gain) and enteric coated (10% theoretical weight gain) according to recommended coating conditions (Colorcon, Table 1).



Table 1. Seal Coating and Delayed Release Coating

O'Hara Labcoat I	Opadry YS-1-7027	Acryl-EZE 93A18597
Tablet Charge (kg)	1	1
Inlet Air Temp. (°C)	70-74	39-42
Drying Air Volume (cfm)	180	125
Tablet Bed Temp. (°C)	43-46	28-32
Exhaust Air Temp. (°C)	52-55	28-32
Atomization Air (psi)	30	9
Pattern Air (psi)	30	15
Spray Rate (g/min.)	15	10
Pan Speed (rpm)	18	20

Table 2 shows the plasticizer type and level (w/w, respect to polymer) and the solids content of the final coating dispersions.

Table 2. Plasticizer Type and Level

Plasticizers (water solubility, 20°C)	Level (% w/w, respect to polymer)	Solids (%)
PEG 8000 (1 in 25)	8, 9, 12, 15	21-22
TEC (1 in 15)	12, 15, 20, 25	22-23
Triacetin (1 in 14)	12, 15, 18, 20, 25, 30	22-24

Sample Analysis

Disintegration and acid uptake testing was performed in a Vankel 35-1200 disintegration apparatus. The disintegration time of the uncoated and coated cores (n=6) in 900ml of 0.1N HCl (pH 1.0), acetate buffer (pH 4.5) and phosphate buffer (pH 6.8) was recorded. The amount of media (HCl acid or acetate buffer) taken up by the coated placebo tablets (acid uptake) was determined by calculating the percent difference between tablet weights before and after exposure to media for 2 hours.


Results and Discussion

Disintegration Results

The uncoated and seal coated cores disintegrated in the acid and buffer phases in less than 1 minute, as expected. The enteric coated placebo tablets remained intact for 2 hours in the acid phases (pH 1.0 and 4.5) and disintegrated in phosphate buffer (pH 6.8) in about 2 minutes, as desired.


Historical data has shown that acid uptake less than 10% corresponds to acceptable enteric protection (no drug release or degradation in the acid phase of dissolution testing). Figure 1 shows that Acryl-EZE 93A plasticized with TEC, PEG 8000 or triacetin had less than 4% acid uptake when tested in 0.1N HCl acid.

Figure 1. Acid Uptake in 0.1N HCl (pH 1.0)

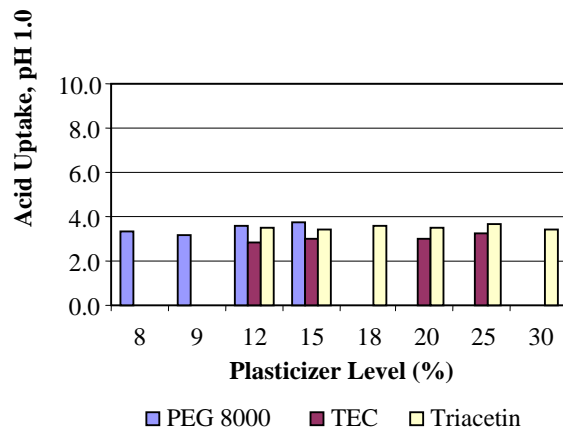

In acetate buffer (pH 4.5), Acryl-EZE 93A plasticized with PEG 8000 had the least acid uptake, followed by TEC and triacetin (Figure 2). PEG 8000 showed a fairly constant acid uptake (about 5%) at each plasticizer level (8-15%). In contrast, TEC and triacetin showed a decrease in acid uptake with an increase in plasticizer level. Specifically, acid uptake decreased from 7.7 to 4.1% as TEC increased from 12 to 25% and it decreased from 10.5 to 5.7% as triacetin increased from 12 to 30%.

Figure 2. Acid Uptake in Acetate Buffer (pH 4.5)

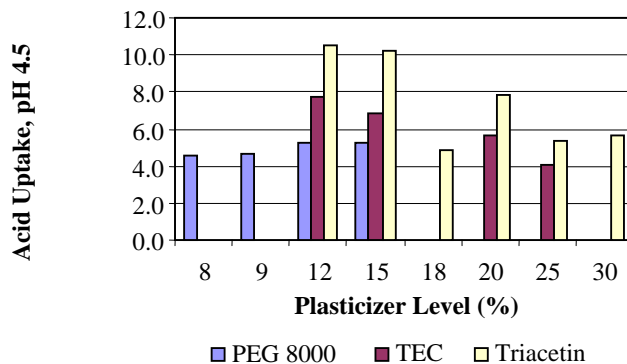

Conclusion

The degree of acid resistance was influenced by the plasticizer type and its concentration in the coating dispersion. All three plasticizers (PEG 8000, TEC, and triacetin) at all tested concentrations within the Acryl-EZE film coating system provided the tablets with enteric protection.

Acryl-EZE 93A provides a flexible pigmented formulation for scientists, allowing them to select the type and level of plasticizer used, while maintaining excellent enteric performance in intermediate pH media.

References

1. Pharmaceutical Coatings Bulletin 102-2, Morflex, Inc., (1994).
2. Felton, L., et al., International Journal of Pharmaceutics 113 (1995) 17-24.
3. Handbook of Pharmaceutical Excipients, Third Edition, American Pharmaceutical Association and Pharmaceutical Press, 2000.
4. Miner, P., et al., American Journal of Gastroenterology 98 (2003) 2616-2620
5. Rohss, K., et al., European Journal of Clinical Pharmacology 60 (2004) 531-539

World Headquarters

Colorcon

415 Moyer Blvd., P.O. Box 24, West Point, PA 19486-0024

Tel: 215-699-7733 Fax: 215-661-2605 Web Site @<http://www.colorcon.com> E.mail: modified_release@colorcon.com

Locations	Telephone	Facsimile	Locations	Telephone	Facsimile
<i>United States</i>			<i>Asia/Pacific</i>		
Santa Ana, California	714-549-0631	714-549-4921	Singapore	65-6438-0318	65-6438-0178
Indianapolis, Indiana	317-545-6211	317-545-6218	Nishiyama, Japan	81-5-4465-2711	81-5-4465-2730
Humacao, Puerto Rico	787-852-3815	787-852-0030	Shanghai, China	86-21-5442-2222	86-21-5442-2229
Quebec, Canada	514-337-8341	514-337-9159	Goa, India	91-832-288-3434	91-832-288-3440
			Seoul, Korea	82-2-2057-2713	82-2-2057-2179
<i>Europe</i>			<i>Latin America</i>		
Dartford, Kent, England	44-1322-293000	44-1322-627200	Buenos Aires, Argentina	54-11-4552-1565	54-11-4552-3997
Massy, France	33-1-6447-9750	33-1-6932-5983	Cotia, Brasil	55-11-4612-4262	55-11-4612-3307
Idstein, Germany	49-6126- 9961- 0	49- 6126-9961-11	Bogota, Colombia	571-418-1202	571-418-1257
Gallarate, Italy	39-0331-776932	39-0331-776831	Santa Fe, Mexico	525-5-3000-5700	525-5-3000-5701
Budapest, Hungary	36-1-200-8000	36 -1-200-8010	Caracas, Venezuela	58-212-442-4819	58-212-442-8724
Barcelona, Spain	34-9-3589-3756	34-9-3589-3792			
Istanbul, Turkey	90-216-465-0360	90-216-465-0361			

The information contained herein, to the best of our knowledge is true and accurate. Any recommendations or suggestions are made without warranty or guarantee, since the conditions of use are beyond our control. Any information contained herein is intended as a recommendation for use of our products so as not to infringe on any patent.

© Colorcon, 2008. The information contained in this document is proprietary to Colorcon and may not be used or disseminated inappropriately. Acryl-EZE[®] and Opadry[®] are registered trademarks of BPSI Holdings LLC.
mr/crs_2007/plasticizer_cy/03.08_rev1