



**SURETERIC®**  
AQUEOUS ENTERIC COATING SYSTEM

Technical Data Sheet  
Coating Application

**Coating Tablets with Sureteric**

Sureteric is a specially blended combination of PVAP (polyvinyl acetate phthalate), plasticizers, and other processing ingredients, specifically designed to meet the enteric coating needs of solid oral dosage forms in the pharmaceutical industry

Coating Parameter	24" Accela-Cota	48" Accela-Cota	60" Accela-Cota
Solvent	Distilled Water	Distilled Water	Distilled Water
Solids content (%w/w)			
<b>Opadry®</b> subcoat	15	15	15
<b>Sureteric®</b>	15	15	15
<b>Opadry II topcoat</b>	15-20	15-20	15-20
Theoretical weight gain (%)			
<b>Opadry subcoat</b>	2	2	2
<b>Sureteric</b>	10	10	10
<b>Opadry II topcoat</b>	3	3	3
Tablet charge (kg) <sup>(1)</sup>	11-13	120-150	280-320
Inlet air temperature (°C)	75-84	70-80	70-80
Drying air volume (m <sup>3</sup> /hr) <sup>(3)</sup>	425+/-42	3227+/-170	4250+/-340
Exhaust air temperature (°C)			
<b>Sureteric</b>	38-41	40-45	40-45
<b>Sub and topcoat</b>	42-44	45-48	45-48
Pre-warm tablet bed (°C)	45-48	45-48	45-48
Spray equipment	2 x Binks 605	4 x Binks 605	5 x Binks 605
Fluid nozzle (mm)	1.8 mm, Binks 66SH	1.8 mm, Binks 66SH	1.8 mm, Binks 66SH
Air cap	External Atom., Oval	External Atom., Oval	External Atom., Oval
Atomizing air pressure (bar) <sup>(3)</sup>	2.4-2.7	4.0-5.5	4.0-5.5
Pattern air pressure (bar) <sup>(3)</sup>	N/A	N/A	N/A
Gun-to-bed distance (cm)	13-18	25-31	25-31
Spray rate (g/min)			
<b>Sureteric</b>	60-70	400-500	700-800
<b>Sub and topcoat</b>	55-60	350-450	650-750
Pan speed (rpm)	14	4-8	3-6
Coating Time (Hours)			
<b>Opadry subcoat</b>	0.5	1	1
<b>Sureteric</b>	2	3-4	3-4
<b>Opadry II topcoat</b>	1	1-1.5	1-1.5

- 1) Maximize pan charge
- 2) Air volume may be decreased to prevent edge chipping
- 3) Adjust to maximize efficiency



### **Additional Considerations**

1. Always maintain a negative air pressure in the pan (more air out than in).
2. Rotate (jog) the table and bed minimally prior to constant rotation, during the spray cycle.
3. After start-up, allow a minimum of 15 minutes for exhaust temperature to equilibrate before making changes in fluid and/or air flows.
4. To maximize enteric performance and adhesion between the sub-coat and enteric interface, reduce the Sureteric spray rate by 15%, for the first 1% weight gain, if any tackiness or sticking is noticed. The delivery rate may then be increased as the cores become sealed with Sureteric.
5. Once Sureteric delivery has begun, maintain a constant flow. Gun needles should remain in an open position for the duration of the Sureteric application.
6. Immediately after applying the topcoat, tablets may be polished using powdered Carnauba wax. Apply no more than 10.0 g per 100 kg of tablets. Apply the wax without exhaust or drying air on the tablets. Run for 5 minutes at 3 rpm, then cool to room temperature with exhaust at low speed for no longer than 5 minutes.

The above parameters are recommended based on Colorcon trial data. Individual product and machine functions may create a need to alter the conditions. Please consult Colorcon if further technical advice is required.

**World Headquarters**

Colorcon

415 Moyer Blvd., P.O. Box 24, West Point, PA 19486-0024

Tel: 215-699-7733 Fax: 215-661-2605 Web Site @<http://www.colorcon.com> E.mail: [modified\\_release@colorcon.com](mailto:modified_release@colorcon.com)

<b>Locations</b>	<b>Telephone</b>	<b>Facsimile</b>	<b>Locations</b>	<b>Telephone</b>	<b>Facsimile</b>
<i>United States</i>			<i>Asia/Pacific</i>		
Santa Ana, California	714-549-0631	714-549-4921	Singapore	65-6438-0318	65-6438-0178
Indianapolis, Indiana	317-545-6211	317-545-6218	Nishiyama, Japan	81-5-4465-2711	81-5-4465-2730
Humacao, Puerto Rico	787-852-3815	787-852-0030	Shanghai, China	86-21-5442-2222	86-21-5442-2229
<i>Europe</i>			Goa, India	91-832-288-3434	91-832-288-3440
Dartford, Kent, England	44-1322-293000	44-1322-627200	Seoul, Korea	82-2-2057-2713	82-2-2057-2179
Bougival, France	33-1-3082-1582	33-1-3082-7879	<i>Latin America</i>		
Idstein, Germany	49-6126-9961-0	49-6126-9961-11	Buenos Aires, Argentina	54-911-4552-1565	54-911-4552-3997
Gallarate, Italy	39-0331-776932	39-0331-776831	Cotia, Brasil	55-11-4612-4262	55-11-4612-3307
Budapest, Hungary	36-1-200-8000	36-1-200-8010	Bogota, Colombia	571-418-1202	571-418-1257
Barcelona, Spain	34-9-3589-3756	34-9-3589-3792	Santa Fe, Mexico	525-5-3000-5700	525-5-3000-5701
Istanbul, Turkey	90-216-465-0360	90-216-465-0361	Caracas, Venezuela	58-212-442-4819	58-212-442-8724

The information contained herein, to the best of our knowledge is true and accurate. Any recommendations or suggestions are made without warranty or guarantee, since the conditions of use are beyond our control. Any information contained herein is intended as a recommendation for use of our products so as not to infringe on any patent.

© Colorcon, 2006. The information contained in this document is proprietary to Colorcon and may not be used or disseminated inappropriately. Sureteric® and Opadry® are registered trademarks of BPSI Holdings Inc. [mr/sureteric/coating\\_with\\_sureterice/10.2006](http://mr/sureteric/coating_with_sureterice/10.2006), Rev 3

First published 07.2003. Revised 10.2006