

## Coating Parameters – Ethanol/Water Solvent Solution

### FOR FORMULATIONS BASED ON HYPROMELLOSE PHTHALATE (HPMC-P)

Opadry® Enteric is a fully formulated, delayed release coating system for pharmaceutical solid oral dosage forms applied by hydro-alcoholic processing techniques. The coating parameters listed in the table below are based on Colorcon data for Opadry Enteric (920 Series) formulations using an ethanol/water solvent solution\*\*. Individual product and machine functions should be taken into account and process parameters adjusted as required.

Opadry Enteric should be applied over tablets seal-coated with a 2%WG of Opadry® (03K19229).

Pan Type		Fully Perforated Side Vented Coating Pan		
Pan Diameter	cm	38.1	60.96	122
	in	15	24	48
Solution Preparation				
Solvent		** ETOH / H2O		
Solids Content	% w/w	12.5	12.5	12.5
Run Objectives				
Theoretical Weight Gain	% w/w	8 – 12	8 – 12	8 – 12
Tablet Charge	kg	1.5 - 3.0	12 – 15	120 – 160
Gun Configuration				
Example Spray Gun	Type	Schlick 931/7-1 S22	Schlick 931/7-1 S35 (ABC)	Schlick 931/7-1 S35 (ABC)
Number of Guns	guns	1	1 – 2	3 – 4
Nozzle Diameter	mm	1.2	1.2	1.2
	in	0.05	0.05	0.05
Gun to Bed Distance	cm	10	17	25
	in	4	7	10
Gun to Gun Distance	cm	N/A	10	15
	in	N/A	4	6
Atomizing Air Pressure	bar	1.4	1.4	1.7
	psi	20	20	25
Pattern Air Pressure	bar	1.4	1.4	1.7
	psi	20	20	25
Process Parameters				
Drying Air Volume	cu.m/hr	250 – 290	424 – 470	2300 – 3060
	cu.ft/min	147 – 175	250 – 275	1350 – 1800
Pan Speed	rpm*	18	14	7
Pre-warm Tablet Bed	°C	33 – 36	33 – 36	33 – 36
	°F	91.4 – 96.8	91.4 – 96.8	91.4 – 96.8
Inlet Air Temperature	°C	36	36	36
	°F	96.8	96.8	96.8
Tablet Bed Temperature	°C	31 – 33	31 – 33	31 – 33
	°F	87.8 – 91.4	87.8 – 91.4	87.8 – 91.4
Spray Rate	g/min	22 – 25	48 – 52	250 – 340***
Approximate Coating Times	min	44 – 195	195 – 338	374 – 550***
*Pan speed should be set so that even mixing of the bed occurs. This will be influenced by tablet size, shape, friability and baffle type in the pan. ** Ethanol/Water mass ratio 80/20*** Estimated				

---

The information contained herein, to the best of Colorcon, Inc.'s knowledge is true and accurate. Any recommendations or suggestions of Colorcon, Inc. with regard to the products provided by Colorcon, Inc. are made without warranty, either implied or expressed, because of the variations in methods, conditions and equipment which may be used in commercially processing the products, and no such warranties are made for the suitability of the products for any applications that you may have disclosed. Colorcon, Inc. shall not be liable for loss of profit or for incidental, special or consequential loss or damages.

Colorcon, Inc. makes no warranty, either expressed or implied, that the use of the products provided by Colorcon, Inc., will not infringe any trademark, trade name, copyright, patent or other rights held by any third person or entity when used in the customer's application.

For more information, contact your Colorcon representative or call:

North America	Europe/Middle East/Africa	Latin America	India	China
<b>+1-215-699-7733</b>	<b>+44-(0)-1322-293000</b>	<b>+54-1-5556-7700</b>	<b>+91-832-6727373</b>	<b>+86-21-61982300</b>

You can also visit our website at [www.colorcon.com](http://www.colorcon.com)



© BPSI Holdings LLC, 2019.

The information contained in this document is proprietary to Colorcon and may not be used or disseminated inappropriately.

All trademarks, except where noted, are property of BPSI Holdings, LLC.

pi\_coat\_par\_opadry\_enteric\_91s\_IPAH2O\_V2\_052019

**This document is valid at the time of distribution. Distributed 25-Mar-2023 (UTC)**